



**ZYMO RESEARCH**

*The Beauty of Science is to Make Things Simple*

# INSTRUCTION MANUAL

## **Quick-DNA™ Viral Kit**

Catalog Nos. **D3015 & D3016**

### **Highlights**

- Quick (*15 minute*) recovery of viral DNA from a wide range of sample sources using Zymo-Spin™ Technology.
- Column design allows DNA to be eluted at high concentrations into minimal volumes ( $\geq 6 \mu\text{l}$ ) of water or elution buffer.
- Eluted DNA is suitable for use in PCR, arrays, and other sensitive downstream applications.

### **Contents**

Product Contents & Specifications.....	1
Product Description.....	2
Buffer Preparation.....	3
Protocol.....	3
Ordering Information.....	4
List of DNA-Related Products.....	5



Satisfaction of all Zymo Research products is guaranteed. If you should be dissatisfied with this product please call 1-888-882-9682.

## Product Contents

<b>Quick-DNA™ Viral Kit (Kit Size)</b>	<b>D3015 (50 preps.)</b>	<b>D3016 (200 preps.)</b>	<b>Storage Temperature</b>
<b>Viral DNA Buffer*</b>	50 ml	2 x 100 ml	Room Temp.
<b>DNA Wash Buffer** (Concentrate)</b>	6 ml	24 ml	Room Temp.
<b>DNA Elution Buffer</b>	4 ml	10 ml	Room Temp.
<b>Zymo-Spin™ IC Columns</b>	50	200	Room Temp.
<b>Collection Tubes</b>	50	200	Room Temp.
<b>Instruction Manual</b>	1	1	-

Note - Integrity of kit components is guaranteed for up to one year from date of purchase. Reagents are routinely tested on a lot-to-lot basis to ensure they provide maximal performance and reliability.

\**Recommended:* Add beta-mercaptoethanol to 0.5%(v/v) i.e., 250 µl per 50 ml or 500 µl per 100 ml.

\*\*Add 24 ml (96 ml for D3016) of 100% ethanol to the DNA Wash Buffer concentrate prior to use.

## Specifications

- **Sample Sources** – Whole blood, serum, plasma, tissue, or cells derived from humans, mice, rabbits, etc.
- **DNA Purity** – High quality DNA is eluted with **DNA Elution Buffer** making it perfect for PCR.  $A_{260}/A_{280} > 1.8$
- **DNA Size Limits** – From 100 bp up to 50 kb.
- **DNA Recovery** – Typically, up to 5 µg total DNA is eluted into 6-10 µl **DNA Elution Buffer**. For DNA 75 bp to 10 kb, the recovery is 70-90%. For DNA 11 kb to 50 kb the recovery is 50-70%.
- **Equipment** – microcentrifuge, vortex

Note - ™ Trademarks of Zymo Research Corporation. This product is for research use only and should only be used by trained professionals. It is not intended for use in diagnostic procedures. Some reagents included with this kit are irritants. Wear protective gloves and eye protection. Follow the safety guidelines and rules enacted by your research institution or facility.


## ZYMO RESEARCH CORP.

The **Genomic DNA Clean & Concentrator™ (D4010, D4011)** can be used to isolate large DNAs >200 kb (i.e., genomic, mitochondrial, viral, etc.) from crude cell lysates including those from Proteinase K digestions.

## Product Description

The **Quick-DNA™ Viral Kit** from Zymo Research provides a streamlined method for rapid isolation of high quality viral DNA from a wide range of biological sources. Denaturation of viral particles and recovery of DNA is accomplished using a unique **Viral DNA Buffer** and **Zymo-Spin™ IC Column**, respectively. Reagents in the kit are optimized for the recovery of viral DNA from whole blood, (fresh and stored), tissue, ascites, cultured cells, and liquid samples. DNA can be eluted with either water or the supplied elution buffer. Eluted DNA is suitable for PCR, arrays, and other downstream applications.

For high-throughput purification (96-well, 5 µg DNA/well), use the **Quick-DNA™ Viral 96 Kit (D3017, D3018)**.

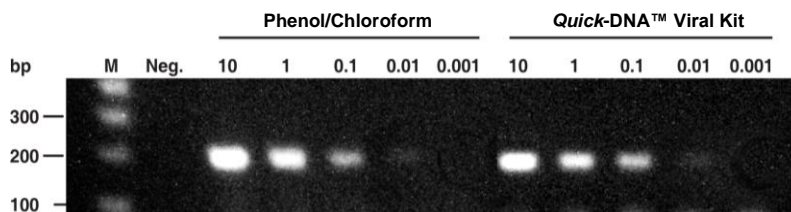

Human HBV DNA was isolated from 10 to 0.001 µl of human serum using phenol/chloroform or Zymo Research's **Quick-DNA™ Viral Kit**. The presence of HBV DNA is evidenced by a ~200 bp PCR product. Lane "M" is a 100 bp DNA Ladder and "Neg" is the negative control for PCR.

For **Technical Assistance**, please contact those at **Zymo Research's Technical Department** at 1-888-882-9682 or E-mail to tech@zymoresearch.com

The **Zymo-Spin™ IC Column** has a maximum capacity of 800 µl. For larger volumes it may be necessary to reload the column.

## **Buffer Preparation**

- ✓ Before starting, add 24 ml 100% ethanol to the 6 ml **DNA Wash Buffer** concentrate (96 ml 100% ethanol to the 24 ml **DNA Wash Buffer** concentrate) to obtain the final **DNA Wash Buffer** solution.
- ✓ *Recommended:* Add beta-mercaptoethanol (user supplied) to the **Viral DNA Buffer** to a final dilution of 0.5%(v/v) i.e., 250 µl per 50 ml or 500 µl per 100 ml.

## **Protocol**

The following protocol is designed for the recovery of viral DNA from ≤200 µl biological liquid (e.g., serum, plasma, CSF, etc.) or ≤200 µl cells in suspension at a concentration of ≤1x10<sup>5</sup> cells/ml. *All centrifugation steps should be performed at 10,000-16,000 x g.*

1. In a 1.5 ml microcentrifuge tube, add 4 volumes of **Viral DNA Buffer** to each volume of sample (e.g., 800 µl **Viral DNA Buffer** to 200 µl serum). Mix briefly by vortexing. Allow to stand at room temperature for 5-10 minutes.

Longer incubation periods may be required for some samples.

2. Transfer the mixture to a **Zymo-Spin™ IC Column** in a **Collection Tube**. Centrifuge for 1 minute. Discard the flow through from the **Collection Tube**.

It may be necessary to centrifuge the sample/Viral DNA Buffer mixture before transferring the supernatant to the Zymo-Spin™ Column to remove particulate matter that may clog the column.

3. Add 300 µl **DNA Wash Buffer** to the column. Centrifuge for 1 minute. Discard the flow through. Repeat this step.

4. Place the **Zymo-Spin™ IC Column** into a microcentrifuge tube. Add 6-10 µl **DNA Elution Buffer** or water directly to the column matrix. Let stand at room temperature for 1 minute. Centrifuge for 1 minute to elute the DNA.

Elution of DNA from the column is dependent on pH and temperature. If water is used, ensure the pH is >6.0. Also, the total yield may be improved by eluting the DNA with higher volumes (20-30 µl) Elution Buffer or water pre-equilibrated to 60-70°C.

DNA is now ready for immediate use or storage.

**Ordering Information**

Product Description	Catalog No.	Kit Size
<b>Quick-DNA™ Viral Kit</b>	D3015	50 preps.
	D3016	200 preps.
<b>Quick-DNA™ Viral 96 Kit</b>	D3017	2 x 96 preps.
	D3018	4 x 96 preps.

For Individual Sale	Catalog No.	Amount
<b>Viral DNA Buffer</b>	D3015-1-50	50 ml
	D3016-1-100	100 ml
<b>DNA Wash Buffer (Concentrate)</b>	D4003-2-6	6 ml
	D4003-2-24	24 ml
<b>DNA Elution Buffer</b>	D3004-4-4	4 ml
	D3004-4-10	10 ml
<b>Zymo-Spin™ IC Columns</b>	C1004-50	50
	C1004-250	250
<b>Collection Tubes</b>	C1001-50	50
	C1001-500	500
	C1001-1000	1,000

**ZYMO RESEARCH CORP.**

## Popular Products From Zymo Research

Product	Description	Kit Size (Preps)	Catalog No. (Format)
<b>Fragment DNA Purification</b>			
<b>DNA Clean &amp; Concentrator™-5</b>	Clean and concentrate up to 5µg DNA into ≥6 µl elution volume in as little as 2 minutes with no wash residue carryover.	50 200 50 200	<b>D4003</b> (uncapped) <b>D4004</b> (uncapped) <b>D4013</b> (capped) <b>D4014</b> (capped)
<b>DNA Clean &amp; Concentrator™-25</b>	Clean & concentrate 25 µg of DNA into ≥25 µl elution volume in as little as 2 minutes with no wash residue carryover.	50 200 50 200	<b>D4005</b> (uncapped) <b>D4006</b> (uncapped) <b>D4033</b> (capped) <b>D4034</b> (capped)
<b>ZR-96 DNA Clean &amp; Concentrator™-5</b>	Quick (15 minute), high-output recovery of up to 5 µg pure DNA into 10-15 µl minimum elution volume allows for highly concentrated DNA.	2x96 4x96	<b>D4023</b> <b>D4024</b>
<b>Genomic DNA Clean &amp; Concentrator™</b>	Quick (5 minute) clean-up of up to 10 µg high molecular weight DNA (≤200 kb) from any enzymatic reaction or impure preparation without precipitations.	25 100	<b>D4010</b> (capped) <b>D4011</b> (capped)
<b>Zymoclean™ Gel DNA Recovery Kit</b>	Purify DNA from high and low-melting agarose gels in minutes	50 200 50 200	<b>D4001</b> (uncapped) <b>D4002</b> (uncapped) <b>D4007</b> (capped) <b>D4008</b> (capped)
<b>ZR-96 Zymoclean™ Gel DNA Recovery Kit</b>	High-throughput DNA purification from high and low-melting agarose gels.	2x96 4x96	<b>D4021</b> <b>D4022</b>
<b>Zymoclean™ Large Fragment DNA Recovery Kit</b>	Purify high molecular weight DNA (≤200 kb) from high and low-melting agarose gels in minutes	25 100	<b>D4045</b> (capped) <b>D4046</b> (capped)
<b>OneStep™ PCR Inhibitor Removal Kit</b>	Fast, one step procedure for removal of PCR inhibitors such as polyphenolics, humic/fulvic acids, melanin, etc. for successful PCR and other downstream applications.	50 2x96	<b>D6030</b> <b>D6035</b>
<b>Plasmid DNA Purification</b>			
<b>Zyppy™ Plasmid Miniprep Kit</b>	Pellet-Free™ plasmid DNA purification in less than 10 minutes. Recover up to 25 µg DNA in as low as 30 µl.	50 100 400	<b>D4036</b> <b>D4019</b> <b>D4020</b>
<b>Zyppy™-96 Plasmid Miniprep</b>	The fastest and simplest high-throughput method for plasmid purification.	2x96 4x96 8x96	<b>D4041</b> <b>D4042</b> <b>D4043</b>
<b>Zyppy™ Plasmid Midiprep Kit</b>	Pellet-Free™ plasmid DNA purification in 15 minutes in a 150 µl minimum elution volume).	25 50	<b>D4025</b> <b>D4026</b>
<b>ZR Plasmid Miniprep™ Classic</b>	Plasmid DNA purification in minutes: (alkaline lysis/spin column format for low 30 µl elution volume).	50 100 400	<b>D4036</b> <b>D4019</b> <b>D4020</b>
<b>Genomic DNA Purification</b>			
<b>Quick-DNA™ Miniprep Kit</b>	Easy purification of genomic DNA from whole blood, plasma, serum, body fluids, buffy coat, lymphocytes, tissue, swabs or cultured cells in as little as 15 minutes <u>without</u> the use of Proteinase K or organic denaturants.	50 200	<b>D3024</b> <b>D3025</b>
<b>Quick-DNA™ 96 Kit</b>	Simple, high throughput (96-well) purification of DNA from whole blood, plasma, serum, body fluids, buffy coat, lymphocytes, tissue, swabs, or cultured cells in about 30 minutes.	2x96 4x96 10x96.	<b>D3010</b> <b>D3011</b> <b>D3012</b>
<b>Environmental DNA Purification Kits</b>	Unique BashingBead™ technology allows isolation of DNA from samples refractory to conventional lysis procedures including tough-to-lyse tissues, soil samples, feces, plants, seeds, insects, bacteria, yeast, filamentous fungi, unicellular and filamentous algae, and protozoa		<b>Visit website for a comprehensive list</b>

Please visit our website to see our complete line-up of products.

**ZYMO RESEARCH CORP.**

Phone: (949) 679-1190 ▪ Toll Free: (888) 882-9682 ▪ Fax: (949) 266-9452 ▪ info@zymoresearch.com ▪ www.zymoresearch.com